



055PR

G-Biosciences ♦ 1-800-628-7730 ♦ 1-314-991-6034 ♦ technical@GBiosciences.com

A Geno Technology, Inc. (USA) brand name

Caspase Substrates & Set

(Cat. #CPS-145, CPS-002, CPS-370, CPS-006, CPS-008,
CPS-009, CPS-013, CPS-113S)



think proteins! think G-Biosciences www.GBiosciences.com

INTRODUCTION

Caspase Substrates & Sets are ready-to-use solutions of caspase substrates for both fluorometric and colorimetric detection of caspase activity in various cell types. These substrates are prepared as 1mM solution and are supplied individually as well as in a set of all caspase substrates for studying multiple apoptosis processes.

ITEMS SUPPLIED

Catalog #	Description	Size
CPS-145	Caspase-1,4,5 Substrate; Ac-WEHD-AFC (1mM)	250µl
CPS-002	Caspase-2 Substrate; Ac-VDVAD-AFC (1mM)	250µl
CPS-370	Caspase-3,7,10 Substrate; Ac-DEVD-AFC (1mM)	250µl
CPS-006	Caspase-6 Substrate; Ac-VEID-AFC (1mM)	250µl
CPS-008	Caspase-8 Substrate; Ac-IETD-AFC (1mM)	250µl
CPS-009	Caspase-9 Substrate; Ac-LEHD-AFC (1mM)	250µl
CPS-013	Caspase-13 Substrate; Ac-LEED-AFC (1mM)	250µl
CPS-113S	Caspase 1-10 and 13 <u>Substrates Set</u>	250µl each

STORAGE CONDITIONS


It is shipped in blue ice. Store at -20°C and is stable for 6-8 months if stored properly.

PREPARATION BEFORE USE

Take out the vial(s) from refrigerator and thaw to RT. Vortex to mix the contents. Avoid repeated freezing and thawing (prepare small aliquots as per requirement). The substrate is ready to use for apoptosis screening and other related assays.

RELATED PRODUCTS

Download our Protein Assays and Bioassays Handbook.

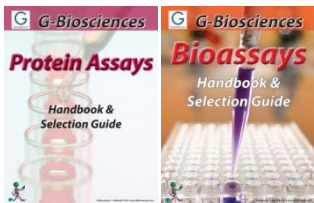

<http://info2.gbiosciences.com/complete-protein-assay-guide>

<http://info2.gbiosciences.com/complete-bioassay-handbook>

For other related products, visit our website at www.GBiosciences.com or contact us.

Last saved: 2/18/2014 TNN



www.GBiosciences.com