IRIS9 - Prestained Protein Ladder



| Catalog Number | Size | Range |
|----------------|--------|---------------|
| PM009-0500 | 500 μΙ | 15 to 180 kDa |

Storage Conditions

Stable for up to 2 weeks at 25°C. Stable for up to 3 months at 4°C. Stable for up to 24 months at -20°C.

Description

The IRIS9 Prestained Protein Ladder is a combination of 9 pre-stained proteins with molecular weights from 15 to 180 kDa. The 9 recombinant proteins are covalently coupled with blue chromophore, while 2 red bands at 70 kDa and 30 kDa, a green band at 15 kDa and a newly designed peacock green band at 60 kDa serve as reference bands. The IRIS9 Prestained Protein Ladder keeps track of the size and separation of proteins during SDS-polyacrylamide gel electrophoresis, approximating the target protein size and validating the Western transfer efficiency on PVDF, nylon, or nitrocellulose membranes.

Kit Content(s)

IRIS9 Prestained Protein Ladder

500µl x 1 vial

Storage Buffer

0.2~0.4 mg/ml of each band of protein is stored in a mixture of solution that contains: 20 mM Trisphosphate at pH 7.5, 2% SDS, 0.2 mM Dithiothreitol, 3.6 M Urea, and 15% (v/v) Glycerol.

Required materials but not provided

- Vertical Electrophoresis system
- Power supplies
- Vortex or equivalent
- Microcentrifuge

Reaction Setup


3μl or 5μl per loading for clear visualization during electrophoresis on 15-well or 10-well mini-gel, respectively. 2.5μl per well for general Western transferring.





Guide for Molecular Weight Estimation (kDa)

Refer to the IRIS9 Prestained Protein Ladder patterns in various electrophoresis conditions:

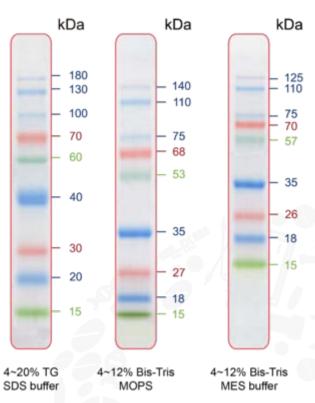

Important notes

The molecular weight of each protein (kDa) was measured against an unstained protein ladder in every electrophoresis condition. Additional data should be considered for a more accurate adjustment.



5F., No. 145, Sec. 3, Beixin Rd., Xindian Dist., New Taipei City, Taiwan (R.O.C.) TEL : +886-2-2462-4956 / FAX : +886-2-2462-8849 www.bio-helix.com